

Newport SR-160-CF

## Fiber Adapted Super Cavity Optical Spectrum Analyzer



Limited Availability  
Used and in Excellent Condition

Open Web Page

<https://www.artisanTG.com/65711-1>

All trademarks, brandnames, and brands appearing herein are the property of their respective owners.



Your **definitive** source  
for quality pre-owned  
equipment.

**Artisan Technology Group**

(217) 352-9330 | [sales@artisanTG.com](mailto:sales@artisanTG.com) | [artisanTG.com](http://artisanTG.com)

- Critical and expedited services
- In stock / Ready-to-ship

- We buy your excess, underutilized, and idle equipment
- Full-service, independent repair center

Artisan Scientific Corporation dba Artisan Technology Group is not an affiliate, representative, or authorized distributor for any manufacturer listed herein.

# Frequency Calibration of the Series 100 and Series 200 SuperCavity™

## Introduction

SuperCavity is a non-confocal, spherically symmetric Fabry Perot interferometer manufactured by Newport Corporation. This compact interferometer has an exceptional efficiency and an extremely high finesse across the specified bandwidth which allows it to be used for detailed investigations and analysis of coherent source mode structures in a near-real time environment.

SuperCavity is not constructed to allow for absolute frequency calibration. However, for many applications, relative calibration of the frequency of the source may be desirable. Before attempting the method of calibration to be described, the incoming beam from the source should be well-aligned and mode-matched to the fundamental cavity mode. (See the Newport Technical Application Note SR-001, "Mode-Matching to the SuperCavity").

A number of methods exist by which the frequency resolving power of Fabry-Perot interferometers such as SuperCavity can be calibrated. The calibration method of Fractional Mode Spacing (FMS) is the subject of this Technical Application Note. Please see the SuperCavity Manual and its references for the Theory of Operation and the theoretical discussion on Fabry-Perot interferometers.

## Basic Equations

The cavity mode structure is composed of both longitudinal and transverse modes that are separated in frequency space (Fig. 1). The mode distribution is given by

$$f_{qmn} = \frac{c}{2L} \left[ q + (m+n+1) \frac{\cos^{-1}(\pm \sqrt{g_1 g_2})}{\pi} \right] \quad (1)$$

where  $c/2L$  = the longitudinal mode spacing,  
 $c$  = the speed of light,  
 $L$  = the cavity length,  
 $m, n$  = transverse mode order = 0,  $\pm 1$ ,  $\pm 2$ , ...,  
 $q$  = longitudinal mode number,  
 $g_i = 1 - L/R_i$ ,  $i = 1, 2$   
 $R_i$  = radius of curvature of the  $i^{\text{th}}$  mirror.

For the Newport SuperCavity, both the front and rear mirrors have radii of 300 mm. The nominal cavity length is either 25.4 millimeters or 20 microns. From Eq. (1) and for  $g_1 = g_2 = g$ , the longitudinal mode spacing is  $c/2L$  and the transverse mode spacing is

$$\text{transverse mode spacing} = \frac{c}{2L} \left( \frac{\cos^{-1} g}{\pi} \right) \quad (2)$$

Eq. (2) may be approximated for  $R \gg L$  by

$$\text{transverse mode spacing} = \frac{c}{\pi \sqrt{2LR}}$$

The ratio of the transverse to the longitudinal mode spacing is referred to as the fractional mode splitting (FMS). The FMS values used to build standard Newport cavities are:

Type of Cavity	Fractional Mode Splitting
SR-100 Series (1")	0.131
SR-200 Series (20μm)	0.00365

With these values and Eq. (2), the length of the respective cavity may be found and used to calculate the longitudinal mode spacing:

Type of Cavity	Longitudinal Mode Splitting $\left( \frac{c}{2L} \right)$
SR-100 Series	$5.984 \pm 0.299$ GHz
SR-200 Series	$7601 \pm 380$ GHz

The inaccuracy in the above mode spacing is due to the corresponding inaccuracy in the cavity length. In absolute frequency space, the transverse mode splitting translates to the following:

Type of Cavity	Transverse Mode Splitting
SR-100 series	$784 \pm 40$ MHz
SR-200 series	$27.7 \pm 1.4$ GHz

The values above refer to the frequency difference between any two transverse modes, and may be used as a means for calibrating the observed spectra. This FMS may further be calibrated by injecting sidebands into the spectrum using an acousto-optic modulator as shown in the SuperCavity Instruction Manual.

For custom FSR SuperCavities, equations (1) and (2) may be used to calculate the FMS values using the appropriate cavity length ( $L$ ).

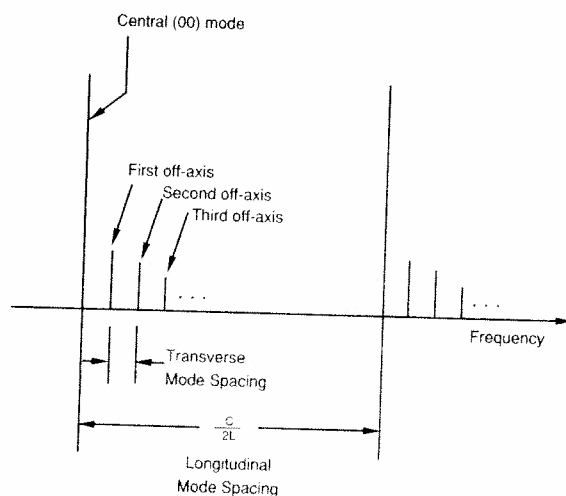


Fig. 1. SuperCavity mode structure and frequency space

DS-11881 (1/12/90)

**U.S. Headquarters:**  
Newport Corporation  
P.O. Box 8020 / 18235 Mt. Baldy Circle  
Fountain Valley, CA 92728-8020  
Telephone: 714-963-9811  
Facsimile: 714-963-2015

**European Headquarters:**  
Federal Republic of Germany  
Newport GmbH  
Telephone: 06151-26116  
Facsimile: 06151-22639

**Canada:**  
Newport Instruments Canada  
Telephone: 416-567-0390  
Facsimile: 416-567-0392

**United Kingdom:**  
Newport Ltd  
Telephone: 0582-769995  
Facsimile: 0582-762655

**Switzerland:**  
Newport Instruments AG  
Telephone: 01-740-2283  
Facsimile: 01-740-2503



**Newport®**

# Artisan Technology Group is an independent supplier of quality pre-owned equipment

## Gold-standard solutions

Extend the life of your critical industrial, commercial, and military systems with our superior service and support.

## We buy equipment

Planning to upgrade your current equipment? Have surplus equipment taking up shelf space? We'll give it a new home.

## Learn more!

Visit us at [artisan<sup>tg</sup>.com](https://www.artisantg.com) for more info on price quotes, drivers, technical specifications, manuals, and documentation.

Artisan Scientific Corporation dba Artisan Technology Group is not an affiliate, representative, or authorized distributor for any manufacturer listed herein.

**We're here to make your life easier. How can we help you today?**

(217) 352-9330 | [sales@artisan<sup>tg</sup>.com](mailto:sales@artisan<sup>tg</sup>.com) | [artisan<sup>tg</sup>.com](https://www.artisantg.com)

