

## Forma Scientific 500mL Erlenmeyer Flask Clamp 4 Inch with Spring and Platform Screw



Limited Availability  
New From Surplus Stock

Open Web Page

<https://www.artisanTG.com/81116-24>

All trademarks, brandnames, and brands appearing herein are the property of their respective owners.



Your **definitive** source  
for quality pre-owned  
equipment.

**Artisan Technology Group**

(217) 352-9330 | [sales@artisanTG.com](mailto:sales@artisanTG.com) | [artisanTG.com](http://artisanTG.com)

- Critical and expedited services
- In stock / Ready-to-ship
- We buy your excess, underutilized, and idle equipment
- Full-service, independent repair center

Artisan Scientific Corporation dba Artisan Technology Group is not an affiliate, representative, or authorized distributor for any manufacturer listed herein.

## Forma Shaker Flask Clip Kits:

Kit Number	Size	Spring Tubes	Screws
236010	25 ml	1	1
236011	50 ml	1	1
236012	125 ml	1	1
236013	250 ml/ 300 ml	1 (w/1 lg. pad)	1
236014	500 ml	1	1
236015	1 L	1	5
236016	2 L	2 (w/10 tubes)	5
236017	4 L	2 (w/10 tubes)	5
236018	6 L	2 (w/12 tubes)	5
236019	2.8 L	2 (w/10 tubes)	5
236028	5 L	2 (w/12 tubes)	5

## Installation

1. Attach the Spring to the Flask Clip Legs as shown in the illustrations.

2. The 2-, 2.8-, 4-, 5-, and 6-Liter Flask Clips use two Springs.

The rubber Spring Tubes are positioned between the Clip Legs as shown in the illustration.


The second Spring is installed around the base of the Flask Clip assembly after it has been mounted to the platform.

3. The cushion pad supplied with the 250 ml/300 ml kit is attached to the topside of the flask clip by peel-off adhesive. The base of the flask rests on the cushion pad during use. (Refer to the illustration below.)


4. Attach the Clip Assembly to the platform with the screws provided.



**Forma Scientific, Inc.**  
**Marietta, Ohio**

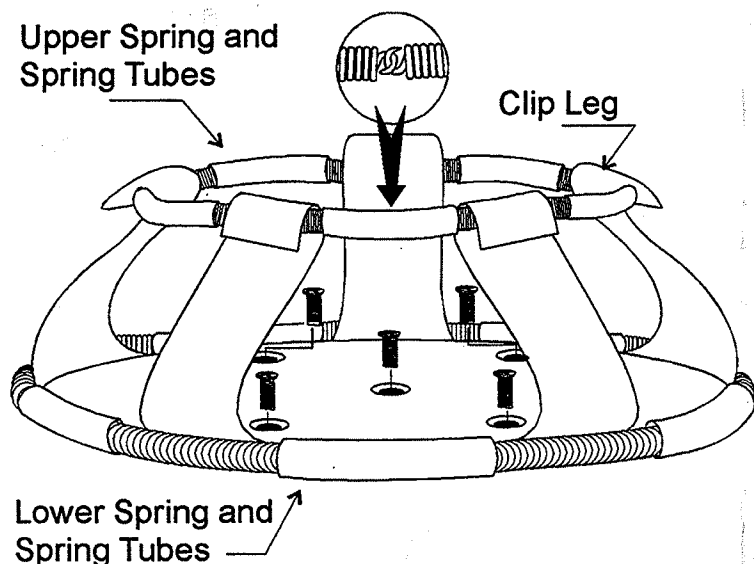

## Flask Clip Kit

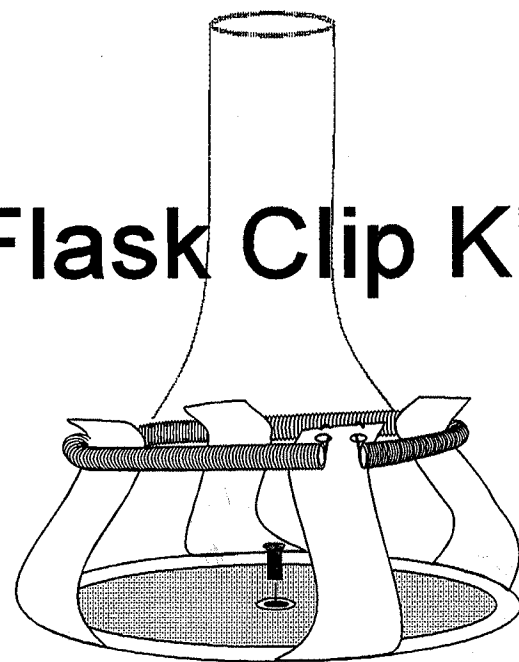

**Instruction Sheet**  
**No. 7236013 Rev. 0**

# Artisan Technology Group is an independent supplier of quality pre-owned equipment

## Gold-standard solutions

Extend the life of your critical industrial, commercial, and military systems with our superior service and support.

## We buy equipment

Planning to upgrade your current equipment? Have surplus equipment taking up shelf space? We'll give it a new home.

## Learn more!

Visit us at [artisan<sup>tg</sup>.com](https://www.artisantg.com) for more info on price quotes, drivers, technical specifications, manuals, and documentation.

Artisan Scientific Corporation dba Artisan Technology Group is not an affiliate, representative, or authorized distributor for any manufacturer listed herein.

**We're here to make your life easier. How can we help you today?**

(217) 352-9330 | [sales@artisan<sup>tg</sup>.com](mailto:sales@artisan<sup>tg</sup>.com) | [artisan<sup>tg</sup>.com](https://www.artisantg.com)

