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CE-DECLARATION OF CONFORMITY
GB
 We declare under our sole responsibility that this product corresponds to the regulations 89/338EEC and 73/232EEC and conforms with the standards or standardized documents EN 61010-1; EN 50081; EN 50082; EN 55011 and EN 60555.

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IKA WERKE GmbH & Co. KG

Staufen, Februar 2004

Wolfgang Buchmann
 Corporate Director - Quality Assurance

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Lasten.

Garantie

You have purchased an original IKA laboratory machine which meets the highest engineering and quality standards.

In accordance with IKA guarantee conditions, the guarantee period is 24 months. For claims under the guarantee please contact

your local dealer. You may also send the machine direct to our

works, enclosing the delivery invoice and giving reasons for the

claim. You will be liable for freight costs.

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4

Contents

Guarantee	2
Safety instructions	4
Correct use	5
Unpacking	5
Commissioning	5
Maintenance and cleaning	5
Technical data	5

The instrument may not be operated in rooms with explosion hazards. Furthermore, it has to be placed upon a fire-proof and/or not burnable surface.

Attention - Magnetism! Effects of the magnetic field have to be taken into account (e.g. data carriers cardiac pavenmakers...).

When using PTFE-coated magnetic bars, the following has to be noted: *Chemical reactions of PTFE occur in contact with molten or dissolved alkaline and alkaline earth metals, as well as with fine-particled powders of metals of the 2. and 3. group of the periodical systems a temperatures above 300-400°C. Only elementary metals do attack it; halogen hydrocarbons have a reversibly swelling effect.*

Source: Römpfs Chemie-Lexikon and „Ullmann“ Vol. 19

Attention: Surface should not be in contact with aggressive medium! Do not place under water!

BIGSQUID0600

Correct use

The IKA BigSquid is a magnetic stirrer suitable for mixing thin-ly fluid media up to 1500 ml by using a magnetic bar up to a length of 30mm. The speed can be adjusted between 0 to a 1500 1/min.

Unpacking

Please unpack the equipment carefully and check for any damages. It is important that any damages which may have arisen during transport are ascertained when unpacking. If applicable a fact report must be set immediately (post, rail or forwarder).
The delivery scope covers:
an IKA BigSquid and operating instructions.

Commissioning

Plug the instrument into a socket with a voltage of 100 ... 240 V \pm 10% and a frequency of 50/60Hz.
Put a magnetic bar into a vessel permeable for magnetic field lines, and fill the vessel with the medium to be processed. Please make sure that the rotary knob for speed adjustment is at the left-hand stop, and then put the vessel on the glass hood. Now you can set the requested speed on the rotary knob.

Technical data

1500ml	Max stirring quantity:
0 ... 1500 1/min	speed range:
100 ... 240 V \pm 10 %	voltage:
50/60 Hz	frequency:
2W	power:
100%	ON duration:
2	protection class:
2	contamination level:
II	overvoltage category:
5 ... 40 °C	ambient temperature:
80%	ambient humidity: (Rel.)
IP 54	Protection to DIN 40050:
800 gr	Weight:
180 x 48mm	Dimensions: (ø x H)

Maintenance and cleaning

The IKA BigSquid is maintenance-free. It is subject only to the natural wear and tear of components and their statistical failure rate. Use only cleansing agents which have been approved by IKA to clean IKA devices. To remove use:
Dyes
isopropyl alcohol
Construction materials
water containing tenside / isopropyl alcohol
Cosmetics
water containing tenside / isopropyl alcohol
Foodstuffs
water containing tenside
Fuels
water containing tenside

For materials which are not listed, please request information from IKA. Wear the proper protective gloves during cleaning of the devices. Electrical devices may not be placed in the cleaning agent for the purpose of cleaning. Before using another than the recommended method for cleaning or decontamination, the user must ascertain with the manufacturer that this method does not destroy the instrument.

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IKAWorks, Inc.

LABORATORY TECHNOLOGY
ANALYZING TECHNOLOGY
PROCESSING EQUIPMENT

North America

IKAWorks, Inc.

2635 North Chase Pkwy SE
Wilmington, NC 28405-7419
USA
TEL. +1 800 733-3037
TEL. +1 910 452-7059
FAX +1 910 452-7693
E-mail: usa@ika.net

IKAWorks, (Asia) Sdn Bhd

LABORATORY TECHNOLOGY
ANALYZING TECHNOLOGY
PROCESSING EQUIPMENT

Asia - Australia

IKAWorks (Asia) Sdn Bhd

No. 17 & 19, Jalan PJU 3/50
Sunway Damansara Technology Park
47810 Petaling Jaya
Selangor, Malaysia
TEL. +60 3 7804-3322
FAX +60 3 7804-8940
E-mail: sales@ika.com.my

IKAJapan Y.K.

LABORATORY TECHNOLOGY
ANALYZING TECHNOLOGY
PROCESSING EQUIPMENT

Japan

IKAJapan Y.K.

293-1 Kobayashi-cho
Yamato Koriyama Shi
639-1026 Japan
TEL. +81 74358-4611
FAX +81 74358-4612
E-mail: japan@ika.de

IKAWorks do Brasil Ltda.

LABORATORY TECHNOLOGY
ANALYZING TECHNOLOGY
PROCESSING EQUIPMENT

South America

IKAWorks do Brasil Ltda.

Estrada do Guernguê, 491
Taquara Jacarepagua, RJ
22713-000 Brasil
TEL. +55 21 2435-9600
FAX +55 21 2435-9601
E-mail: brasil@ika.net

IKAWorks Guangzhou

LABORATORY TECHNOLOGY
ANALYZING TECHNOLOGY
PROCESSING EQUIPMENT

China

IKAWorks Guangzhou

173-175 Friendship Road
Guangzhou Economic & Technological
Development Zone
Guangzhou 510730, P.R.CHINA
TEL. +86 20 8222-6772
FAX +86 20 8222-6776
E-mail: sales@ikagz.com.cn

Artisan Technology Group is your source for quality new and certified-used/pre-owned equipment

- FAST SHIPPING AND DELIVERY
- TENS OF THOUSANDS OF IN-STOCK ITEMS
- EQUIPMENT DEMOS
- HUNDREDS OF MANUFACTURERS SUPPORTED
- LEASING/MONTHLY RENTALS
- ITAR CERTIFIED SECURE ASSET SOLUTIONS

SERVICE CENTER REPAIRS

Experienced engineers and technicians on staff at our full-service, in-house repair center

*InstraView*SM REMOTE INSPECTION

Remotely inspect equipment before purchasing with our interactive website at www.instraview.com ↗

WE BUY USED EQUIPMENT

Sell your excess, underutilized, and idle used equipment. We also offer credit for buy-backs and trade-ins. www.artisanng.com/WeBuyEquipment ↗

LOOKING FOR MORE INFORMATION?

Visit us on the web at www.artisanng.com ↗ for more information on price quotations, drivers, technical specifications, manuals, and documentation

Contact us: (888) 88-SOURCE | sales@artisanng.com | www.artisanng.com