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**CE-DECLARATION OF CONFORMITY**

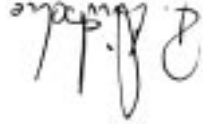
**GB**  
We declare under our sole responsibility that this product corresponds to the regulations 89/36EEC and 73/023EEC and conforms with the standards or standardized documents EN 61 010; EN 50 082; EN 55 014; EN 55 114 and EN 60 555.

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Staufen, Mai 2002

Reiner Diezsch  
Vice President

Wolfgang Buchmann  
Corporate Director - Quality Assurance

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**Garantie**

You have purchased an laboratory machine which meets the highest engineering and quality standards.  
In accordance with guarantee conditions, the guarantee period is 24 months. For claims under the guarantee please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

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2,4 kg

5

The instrument may not be operated in rooms with explosive atmosphere. Effects of the magnetic field have to be taken into account (e.g. data carriers, cardiac pacemakers ...).

When using PTFE-coated magnetic bars, the following have to be noted: Chemical reactions of PTFE occur in contact with molten or dissolved alkaline and alkaline earth metals, as well as with fine-particled powders of metals of the 2. and 3. group of the periodic system at temperatures above 300-400°C.

Only elementary fluorine, chlorine trifluoride and alkaline metals do attack PTFE, halogen hydrocarbons have a reversibly swelling effect.

Source: Römpps Chemie-Lexikon and "Ullmann" Bd.19

### Safety instructions

Guarantee	2
Safety instructions	6
Proper use	6
Unpacking	6
Commissioning	7
Switching on	7
Function Stirring	7
Maintenance and Cleaning	7
Associated standards and regulations	8
Accessories	8
Technical data	8
Information for care and maintenance of the heating plate vers. C	12
List of spare parts REO basic C	14
Spare parts diagram REO basic C	15

### Contents

Even in the case of repair work, the unit must only be opened by a qualified engineer. The mains plug must be pulled before opening the unit. The voltage-carrying components inside the unit may continue to carry a voltage for some time after the mains plug is pulled.

### Proper use

The IKAMAG REO basic C is a laboratory device. It is used in laboratories in the chemical industry, schools and pharmacies to name a few. The devices are suitable for mixing of substances which are placed on mounting plates. The mixing intensity depends on the motor speed and the size of the magnetic stirring bar.

### Unpacking

Please unpack the equipment carefully and check for any damages. It is important that any damages which may have arisen during transport are ascertained when unpacking. It applies cable a fact report must be set immediately (post, rail or forwarder).

The delivery scope covers: an REO C, connecting cable and operating instructions.

REO BC 0502

The REO basic C is maintenance-free. It is subject only to the natural wear and tear of components and their statistical failure rate. **When ordering spare parts, please give the manufacturing number shown on the type plate, the machine type and the name of the spare part.** Please send in equipment for repair only after it has been cleaned and is free from any materials which may constitute a health hazard. Use only cleansing agents which have been approved by IKA to clean IKA devices. To remove use:  
 Dyes isopropyl alcohol  
 Construction materials water containing tenside / isopropyl alcohol  
 Cosmetics water containing tenside / isopropyl alcohol  
 Foodstuffs water containing tenside  
 Fuels water containing tenside  
 For materials which are not listed, please request information from IKA. Wear the proper protective gloves during cleaning of the devices. Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.  
 Before using another than the recommended method for cleaning or decontamination, the user must ascertain with the manufacturer that this method does not destroy the instrument.

## Maintenance and Cleaning

The function stirring is started and stopped by its corresponding switch. If the function "Stirring" is started, the LED "Mot" describes the the imprint "Mot" shines green.  
 The speed of the drive can be adjusted on the "Mot" dial with the corresponding scale from 50 to 1500 rpm.  
 The motor speed of the instrument is kept constant by a control circuit.

## Function Stirring

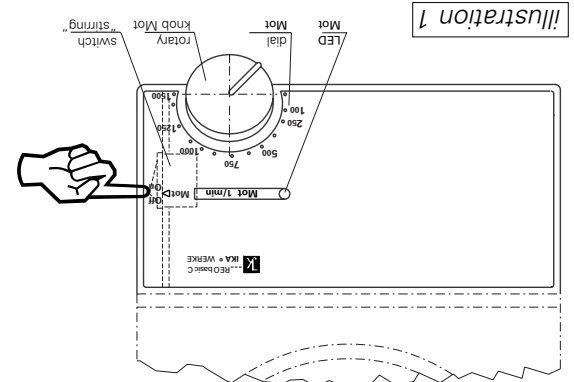
Illustration 1  
 After the instrument has been switched on, it is in "standby"-operation. The function "stirring" is started with the help of switch on the right bottom of the housing. (see illustration 1)

## Switching on

Check whether the voltage specified on the type plate matches the mains voltage available. The power socket used must be earthed (protective earth conductor contact). If these conditions are met, the device is ready to operate after plugging in the mains plug. If these procedures are not followed, safe operation cannot be guaranteed and/or the equipment may be damaged.  
 Observe the ambient conditions (temperature, humidity, etc.) listed under Technical Data.

## Commissioning

Protective conductor.  
On the under side of the mounting plate are areas (bare metal) and can be used as a test point for the protective conductor. It may be necessary to remove an oxide coating for the subsequent tests.

### Associated standards and regulations

**Construction in accordance with the following safety standards**  
EN 61 010-1  
UL 3101-1  
CAN/CSA C22.2 (1010-1)

**Construction in accordance with the following EMC standards**  
EN 61 326-1+A1

**Associated EU guidelines**  
EMC-guidelines: 89/336/EWG  
Machine guidelines: 73/023/EWG

### Accessories

**RS 1** set of stirring bars  
**H16V** support rod  
**RSE** PTFE stirring bar remover  
**H 44** cross sleeve  
**H 36** holding rod

**Recommended IKA-stirrers**  
**TRIKA**-stirrers PTFE, triangular lengths 25, 40 and 55 mm  
**IKAFILON**-stirrers PTFE lengths 10 and 15 mm dia 6mm lengths 20, 25, 30, 40, 50 and 60 mm dia 8mm

### Technical data

design voltage: **VAC** 220 - 240 ±10%  
design frequency: **Hz** 50/60  
order **VAC** 115 ±10%  
order **VAC** 100 ±10%

**Drive:** EC-motor with ball bearings, without brushes  
Max. power consumption **W** 12 / 5  
speed range: / output: **rpm** 50...1500  
infinately variable setting  
speed variety (without load, **%** -5.....+15  
1500rpm):

string quantity (H<sub>2</sub>O): **l** 20  
diameter of mounting plate: **mm** 135  
permitted ambient temperature: **°C** +5 to +40  
permitted relative humidity: **%** 80  
permitted duration of operation: **%** 100  
protection class acc. DIN 40 050: **IP** 42

**fuses:** F1/F2 10AT/250V  
Id-Nr 14 080 00  
contamination level: **2**  
overvoltage category: **II**  
protection class: I(protective earth)

operation at a terrestrial altitude **m** max 2000 above sea level  
dimensions: **mm** 160 x 280 x 85  
weight: **kg** 2,4

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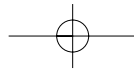
## Information for Care and Maintenance of the Mounting Plate with Technical Enamel Coating





The technical enamel coating makes the mounting plate easier to care for and more resistant to acids and bases. Because of it, however, the mounting plate is also more susceptible to extreme fluctuations in temperature and the force of impact. This can result in cracks forming or the coating flaking off.

Make certain that the bottom of the placing vessel is even, clean and dry. The bottom of the placing vessel must not have any sharp grooves, sides or edges. Remove residues of bases and acids immediately. We recommend most strongly that you clean the mounting plate regularly.

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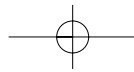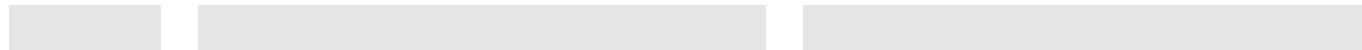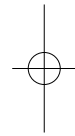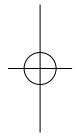14

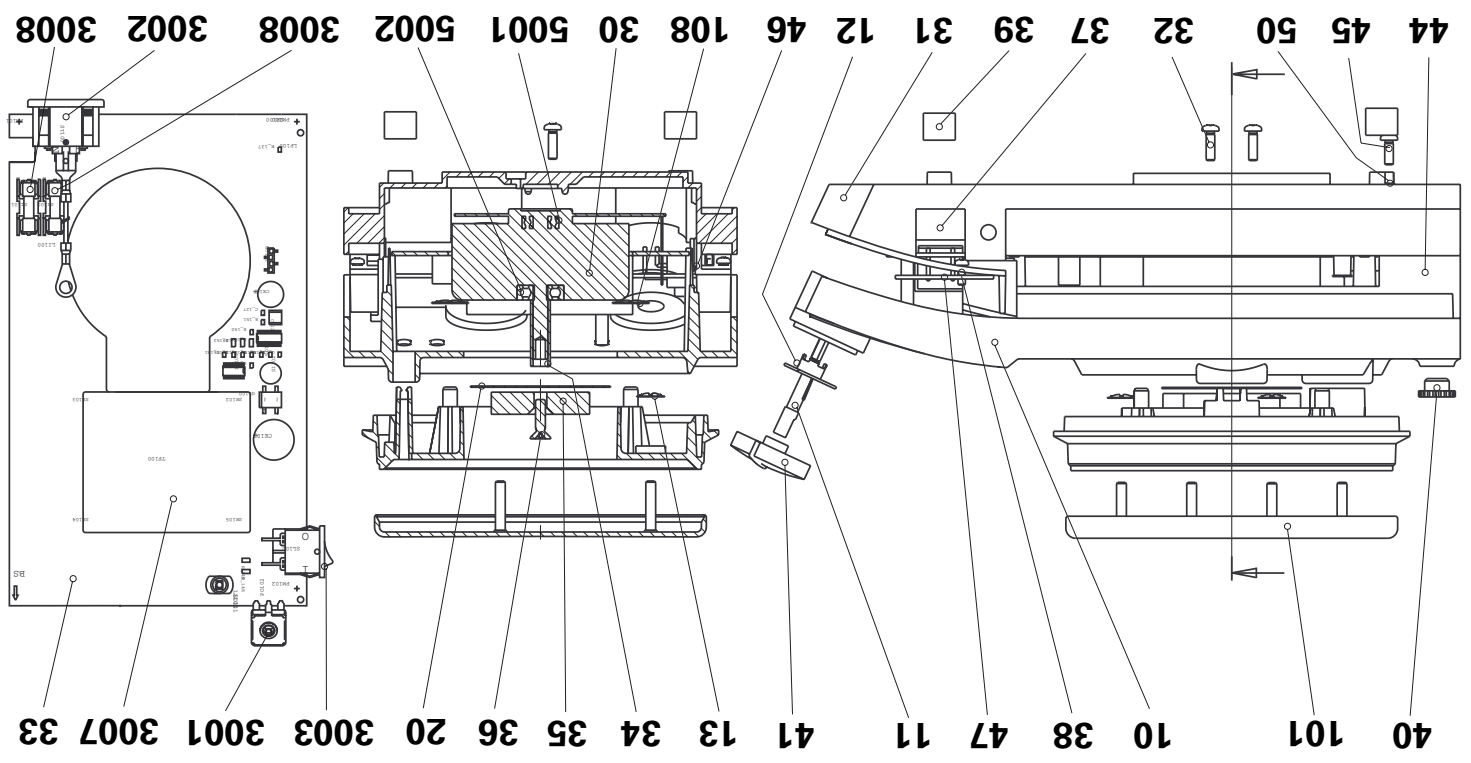
**List of spare parts REO basic C**

Nr.	Designation
10	housing-upper part
11	potentiometer shaft
12	clamping bushing
13	spring nut
20	washer
30	EC - motor
31	housing-bottom part
32	plastic screw
33	PCB control
34	bushing
35	magnet
36	countersunk screw
37	plastic cover left
38	tapfite countersunk screw
39	cap
40	screwed plug
41	operation knob
44	cover
45	tapfite chese-head screw
46	plastic cover right
47	fixing plate
50	lock washer
101	mounting plate
108	spring nut
3001	potl 1k
3002	electric mains
3003	rocker switch motor
3007	trafo
3008	fuse 6,3AT
3010	connector 3pol
3011	connector 4pol
5001	grooved ball bearing
5002	grooved ball bearing

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**IKAWORKS, INC.**

LABORATORY TECHNOLOGY  
ANALYZING TECHNOLOGY  
PROCESSING EQUIPMENT

**IKAWorks, (Asia) Sdn Bhd**

LABORATORY TECHNOLOGY  
ANALYZING TECHNOLOGY  
PROCESSING EQUIPMENT

The Global Market of Laboratory Equipment  
<http://www.labworld-online.com>

Europa - Afrika

**IKAWERKE GMBH & CO.KG**  
Janke & Kunkel-Str. 10  
D 79 219 STAUFEN  
GERMANY  
TEL. 07633/831-0  
FAX 07633/831-98  
E-mail: sales@ika.de  
Internet: <http://www.ika.net>

America

**IKAWORKS, INC.**  
2635 NORTH CHASE PKWY. SE  
WILMINGTON, NC 28405-7419  
TEL. 800/733-3037  
TEL. 910/452-7059  
FAX 910/452-7693  
E-mail: usa@ika.net

Asien - Australien

**IKAWorks (Asia) Sdn Bhd**  
(Company No. 340448-K)  
Lot 2, Jalan Indah 1/2  
Taman Industri Rawang Indah  
48000 Rawang  
Selangor, Malaysia  
TEL: (603) 6093 3322  
FAX: (603) 6093 3940  
E-mail: [ika@tm.net.my](mailto:ika@tm.net.my)

## Artisan Technology Group is your source for quality new and certified-used/pre-owned equipment

- FAST SHIPPING AND DELIVERY
- TENS OF THOUSANDS OF IN-STOCK ITEMS
- EQUIPMENT DEMOS
- HUNDREDS OF MANUFACTURERS SUPPORTED
- LEASING/MONTHLY RENTALS
- ITAR CERTIFIED SECURE ASSET SOLUTIONS

### SERVICE CENTER REPAIRS

Experienced engineers and technicians on staff at our full-service, in-house repair center

### *InstraView*<sup>SM</sup> REMOTE INSPECTION

Remotely inspect equipment before purchasing with our interactive website at [www.instraview.com](http://www.instraview.com) ↗

### WE BUY USED EQUIPMENT

Sell your excess, underutilized, and idle used equipment. We also offer credit for buy-backs and trade-ins. [www.artisanng.com/WeBuyEquipment](http://www.artisanng.com/WeBuyEquipment) ↗

### LOOKING FOR MORE INFORMATION?

Visit us on the web at [www.artisanng.com](http://www.artisanng.com) ↗ for more information on price quotations, drivers, technical specifications, manuals, and documentation

**Contact us:** (888) 88-SOURCE | [sales@artisanng.com](mailto:sales@artisanng.com) | [www.artisanng.com](http://www.artisanng.com)