



## Artisan Technology Group is your source for quality new and certified-used/pre-owned equipment

- FAST SHIPPING AND DELIVERY
- TENS OF THOUSANDS OF IN-STOCK ITEMS
- EQUIPMENT DEMOS
- HUNDREDS OF MANUFACTURERS SUPPORTED
- LEASING/MONTHLY RENTALS
- ITAR CERTIFIED SECURE ASSET SOLUTIONS

### SERVICE CENTER REPAIRS

Experienced engineers and technicians on staff at our full-service, in-house repair center

### *InstraView*<sup>SM</sup> REMOTE INSPECTION

Remotely inspect equipment before purchasing with our interactive website at [www.instraview.com](http://www.instraview.com) ↗

### WE BUY USED EQUIPMENT

Sell your excess, underutilized, and idle used equipment. We also offer credit for buy-backs and trade-ins. [www.artisanng.com/WeBuyEquipment](http://www.artisanng.com/WeBuyEquipment) ↗

### LOOKING FOR MORE INFORMATION?

Visit us on the web at [www.artisanng.com](http://www.artisanng.com) ↗ for more information on price quotations, drivers, technical specifications, manuals, and documentation

**Contact us:** (888) 88-SOURCE | [sales@artisanng.com](mailto:sales@artisanng.com) | [www.artisanng.com](http://www.artisanng.com)

# POLYTRON®

DISPERSING AND MIXING TECHNOLOGY BY KINEMATICA

## Dispersing Technology for the Laboratory

POLYTRON® Systems PT 1200 B/C/CL - Handheld Homogenizers



# KINEMATICA AG

Dispersing and Mixing Technology

# Handy, powerful and safe

## A whole family for small amounts

With the guiding principle „No compromises on small working volumes“, the POLYTRON® PT 1200 handheld disperser family offers three powerful and safe versions. Equipped with different POLYTRON® Dispersing-Aggregates they are suitable for dispersing, homogenizing, emulsifying, suspending, blending, etc. small volumes between approx. 0.1 ml (e.g. Eppendorf tubes) and approx. 250 ml.

### PT 1200 B - independent

- accumulator drive with 30 Watt Leistung
- two eligible speeds 15000 / 25000 rpm
- three exchangeable dispersing aggregates
- quick coupling for dispersing aggregates
- 1h - quick battery charger

### PT 1200 C - safe

- powerful 100 Watt drive unit
- safety transformer with tension outlet 18V
- variable speeds up to approx. 22000 rpm (depending on dispersing aggregate and media)
- three exchangeable dispersing aggregates
- quick coupling for dispersing aggregates
- splash proof housing

### PT 1200 CL - economic


- powerful 100 Watt drive unit
- transformer with tension outlet 18V
- variable speeds up to approx. 22000 rpm (depending on dispersing aggregate and media)
- three exchangeable dispersing aggregates
- quick coupling for dispersing aggregates

### Quick coupling

POLYTRON® handheld dispersers are not only sturdy, safe and powerful but also convincing by their handiness. A quick coupling system allows an aggregate change even with one hand. No screwing, no searching for correct position ...

### Dispersing aggregates in Easy-Clean design

As a standard POLYTRON® dispersing aggregates have a double bearing and are made of finest stainless steel (1.4435/316L). As a surplus the dispersing aggregates for the PT 1200 family are equipped with saw teeth that increases their efficiency in preliminary reduction of the sample and considerably reduces the processing time. The teeth are secured by a safety ring which prevents the teeth from bending. Dispersing aggregates in EC design, EC stands for *EasyClean™*, are equipped with special slide bearings made from high quality plastics which convince with excellent mechanical characteristics and superior chemical resistance. They can be disassembled with effortless ease, are easily cleaned and can be sterilised and autoclaved.

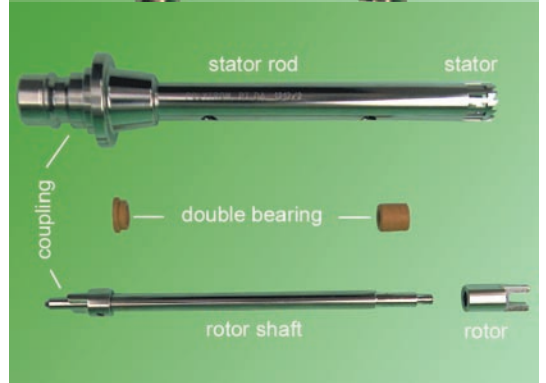

# POLYTRON® PT 1200 family- Overview

Dispersing aggregate / model	PT-DA 1205/2EC	PT-DA 1207/2EC	PT-DA 1212/2EC
<b>Description</b>	suitable for smallest amounts and Eppendorf tubes	suitable for small amounts and test tubes	suitable for small and larger amounts
<b>Shaft length</b>	100 mm	105 mm	105 mm
<b>Stator-Ø</b>	5.5 mm	7.5 mm	12 mm
<b>Rotor-Ø</b>	3 mm	5 mm	9 mm
<b>Tip speed (max.)</b>	3.5 m/s	6 m/s	11 m/s
<b>Working volume</b>	approx. 0.1 - 5 ml	approx. 0.3 - 10 ml	approx. 2 - 250 ml
<b>Recommended dispersing vessel</b>	KINEMATICA dispersing vessels from GS 15 upwards or a vessel with min. 6 mm inside diameter	KINEMATICA dispersing vessels from GS 15 upwards or a vessel with min. 8 mm inside diameter	KINEMATICA dispersing vessels from GS 15 upwards or a vessel with min. 15 mm inside diameter
<b>Applications</b>	<ul style="list-style-type: none"> <li>• Dispersing smallest amounts of plant, human or animal tissues</li> <li>• Dissolving of solids</li> <li>• Extractions</li> </ul>	<ul style="list-style-type: none"> <li>• Dispersing of cells and tissues in small amounts of liquid</li> <li>• Dissolving of solids</li> <li>• Extractions</li> </ul>	<ul style="list-style-type: none"> <li>• Dispersing of plant, human or animal tissues, resins, pigments</li> <li>• Extractions</li> <li>• Producing of suspensions / emulsions</li> </ul>
<b>Order Code</b>	11030005	11030020	11030040



Drive unit / Model	Order Code
PT 1200 B with battery charger 230 V	11010020
PT 1200 B with battery charger 115 V	11010021
Quick battery charger 230 V (approx. 1h)	11080005
Quick battery charger 115 V (approx. 1h)	11080006
PT 1200 C Europe 230 V, 50/60 Hz	11010010
CH 230 V, 50/60 Hz	11010011
Others 115 V, 50/60 Hz	11010012
PT 1200 C 230 V, 50/60 Hz	11010015
115 V, 50/60 Hz	11010016



Accessories		
AP 1200	Zubehörpaket mit Basisstation und Stativ inkl. Transportkoffer	11040005
GS 15	Dispersing vessel for 35 ml	11050010
GS 25	Dispersing vessel for 150 ml	11050011
GS 40	Dispersing vessel for 500 ml	11050012



Other dispersing and homogenizing vessels and accessories on request



# Finest Solutions ... worldwide !

A wide range of products with the finest solution for every application, matched by the superior quality of all components make KINEMATICA a world leader for Dispersing and Mixing technology.

## With Keen Eyes ...

Development knows no rest at KINEMATICA. Our systems are constantly upgraded to reflect the state-of-the-art technology and the needs of the market. KINEMATICA has always had a keen eye for major advances, as proved by various patents. The rotor/stator principle for dispersing technology was conceived and developed at KINEMATICA and has become the world standard.

## ... For Your Needs

Applications for Dispersing/Homogenizing Technology are always on the increase. KINEMATICA continually creates new areas of processing applications based on innovations. We do not neglect the human factor as it is exemplified by the noise reduction of our high-speed systems.

Please, get in touch with us, our local appointed agent or your preferred dealer. We will be pleased to supply more details or arrange for a personal demonstration of our POLYTRON® and MEGATRON® systems.

# Feinste Lösungen ... weltweit !

Eine breite Produktpalette mit der optimalen Lösung für jede Applikation, sowie der hohe Qualitätsstandard aller Systemteile machen KINEMATICA zum weltweit führenden Unternehmen im Bereich der Rotor/Stator-Dispergier- und Mischtechnik.

## Mit offenen Ohren...

KINEMATICA entwickelt die Systeme laufend weiter und bringt sie konzeptionell und konstruktiv auf den neuesten Stand der Technik. Denn KINEMATICA hat die Ohren noch nie vor dem verschlossen, was unerhörte Fortschritte bringt - davon zeugen auch verschiedene Patente.

So hat sich das von KINEMATICA entwickelte Rotor/Stator-Dispergierprinzip weltweit durchgesetzt und bewährt.

## ... für Ihre Bedürfnisse

Die Einsatzbereiche der Rotor/Stator-Dispergiertechnik erweitern sich ständig. Das nicht zuletzt deshalb, weil KINEMATICA durch Innovationen, wie erhöhte Umfangsgeschwindigkeiten, neue Arbeitsgebiete und neue Applikationen erschliesst - ohne dabei menschliche Aspekte zu vernachlässigen. Erinnerung sei an das Bemühen um die Geräuschkürzung am Arbeitsplatz.

Wenden Sie sich an uns. Gerne senden wir Ihnen weitere Unterlagen oder führen Ihnen unsere POLYTRON®- und MEGATRON®-Systeme vor.

Represented by / Vertreten durch:

**POLYTRON**  
**BIOTRONA**  
**MEGATRON**  
**REACTRON**  
**POLYMIX**

Dispersing Technology for Laboratory and Pilot Plant  
Batch- and In-line-Dispersing equipment

**KINEMATICA AG**  
Dispergier- und Mischtechnik

**POLYTRON**  
**BIOTRONA**  
**MEGATRON**  
**REACTRON**  
**POLYMIX**

Homogenizing- and Dispersing Technology  
In-line- und Batch-Processors for Pilot Plant- and Production

**KINEMATICA AG**  
Dispergier- und Mischtechnik



# KINEMATICA AG

Dispersing- and Mixing Technology

Luzernerstrasse 147a  
CH-6014 Littau / Lucerne  
Switzerland

Tel.  
Fax  
e-Mail  
web

+41-41-259 65 65  
+41-41-259 65 75  
laboratory@kinematica.ch  
www.kinematica.ch





## Artisan Technology Group is your source for quality new and certified-used/pre-owned equipment

- FAST SHIPPING AND DELIVERY
- TENS OF THOUSANDS OF IN-STOCK ITEMS
- EQUIPMENT DEMOS
- HUNDREDS OF MANUFACTURERS SUPPORTED
- LEASING/MONTHLY RENTALS
- ITAR CERTIFIED SECURE ASSET SOLUTIONS

### SERVICE CENTER REPAIRS

Experienced engineers and technicians on staff at our full-service, in-house repair center

### *InstraView*<sup>SM</sup> REMOTE INSPECTION

Remotely inspect equipment before purchasing with our interactive website at [www.instraview.com](http://www.instraview.com) ↗

### WE BUY USED EQUIPMENT

Sell your excess, underutilized, and idle used equipment. We also offer credit for buy-backs and trade-ins. [www.artisanng.com/WeBuyEquipment](http://www.artisanng.com/WeBuyEquipment) ↗

### LOOKING FOR MORE INFORMATION?

Visit us on the web at [www.artisanng.com](http://www.artisanng.com) ↗ for more information on price quotations, drivers, technical specifications, manuals, and documentation

**Contact us:** (888) 88-SOURCE | [sales@artisanng.com](mailto:sales@artisanng.com) | [www.artisanng.com](http://www.artisanng.com)